



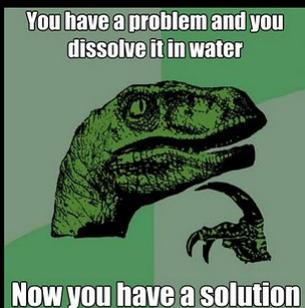

Solutions

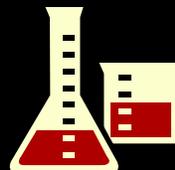

Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

Terms

Solvent

Single substance that does the dissolving
substance present in the largest amount

Solute

1 or more substance that is dissolved
substance present in the lower amount

Solution

The result of dissolving the solute in a solvent

Solubility

Quantity of a solute that will dissolve at a fixed temperature
Typically expressed a grams solute/per 100 (mL or g)



Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

Terms

Saturated

Solution contains the maximum amount of solute
A dynamic equilibrium exists



Unsaturated

Solution contains less than the maximum amount of solute

Supersaturated


Solution contains more than the maximum amount of solute
Carefully prepared
Unstable



Copyright Larry P. Taylor, Ph.D. All Rights Reserved



LPT

Solubility Curve

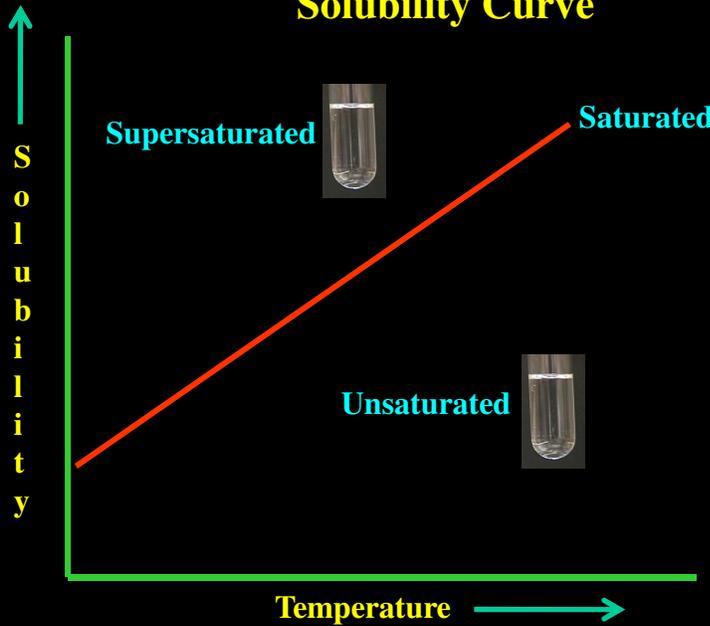


Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

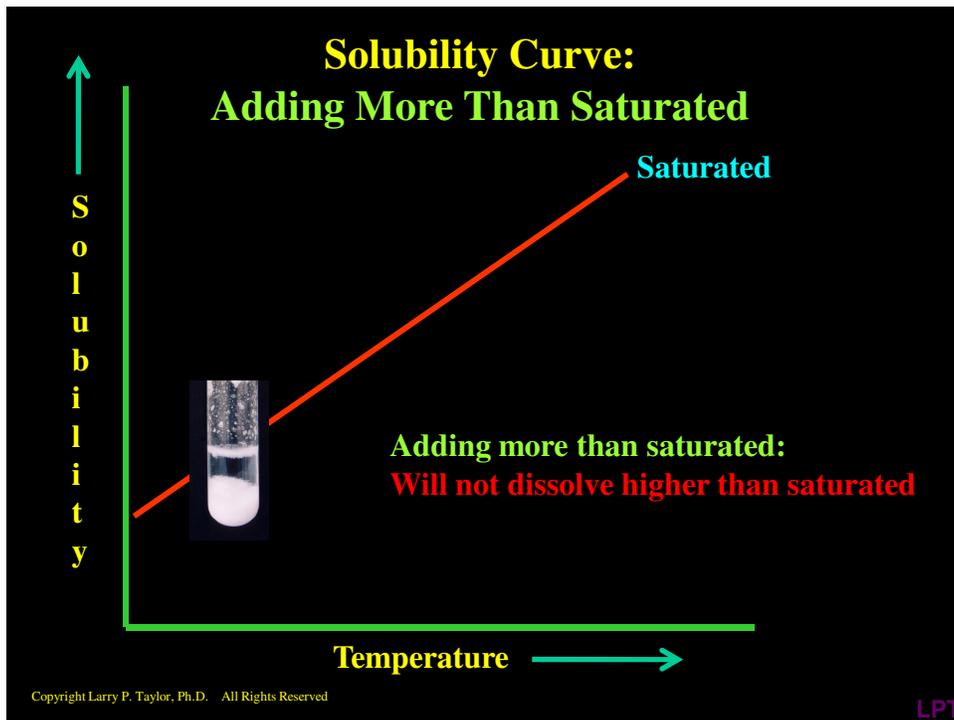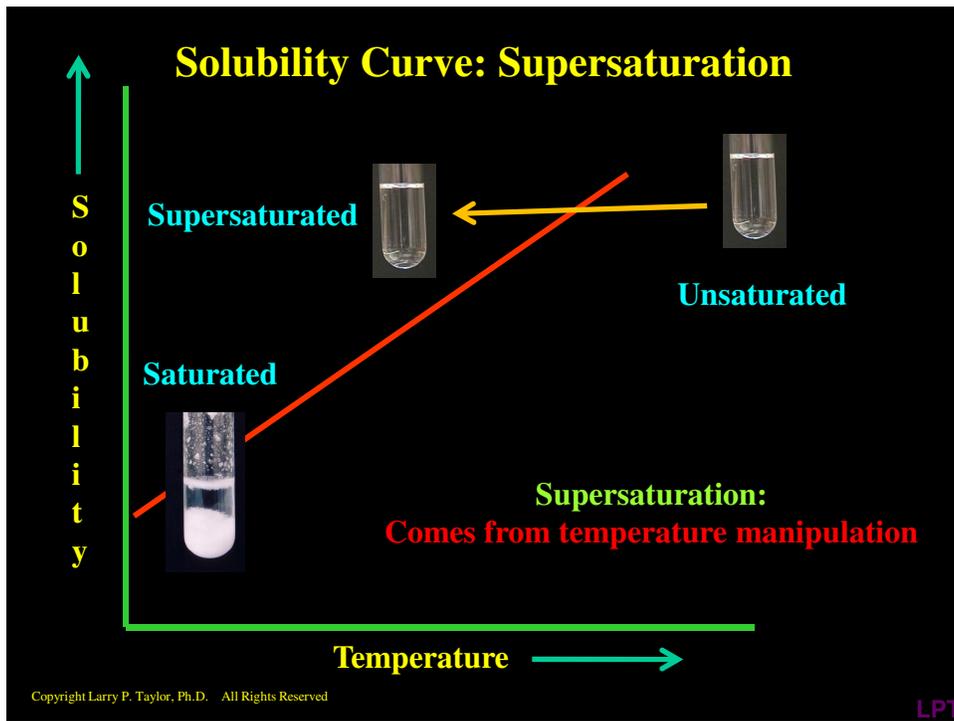

Solubility Curves

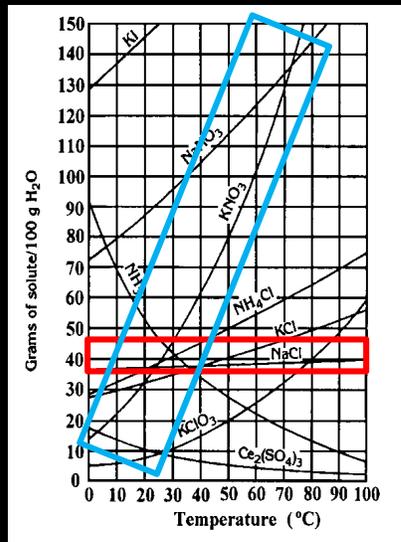

Plot of Solubility vs. Temperature

For each compound (line):

Saturated
On the solubility line


Unsaturated
Below the line

Supersaturated
Above the line

Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

Solubility Curves

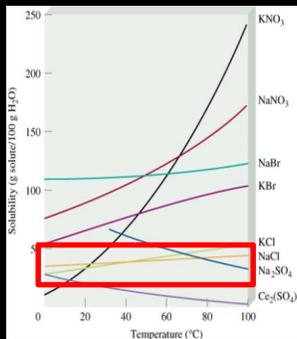

NaCl curve is "flat line"

Solubility does not change with temperature

Consider this:

Our nervous system is dependent on Na⁺



If Na⁺ solubility changed with temperature,
Our nervous systems would alter with temperature

Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT





Solutions Alter Physical Properties Dependent on Concentration of Solute



$$\Delta T = K_b m$$

$$\Delta T = K_f m$$

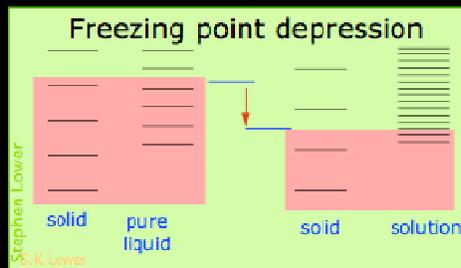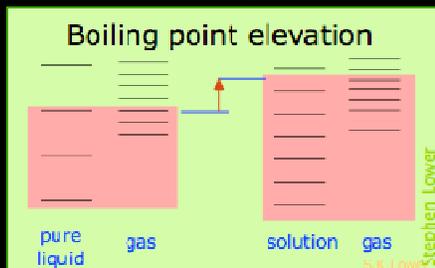

Adding salt to water:

Raises Boiling Point

Raises Cooking Temperature

(but not enough to be significant)

Adding salt to water:

Lowers Freezing Point

Keeps water on sidewalks liquid

Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

WATER ON MARS

It has been found

www.jacanaent.com




Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT



Solutions Lab

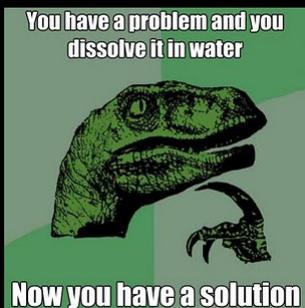

Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

Online Lab




Purpose

Observe the solubility- temperature relationship
Isolate 1.00 gram of NaCl by evaporation

Procedure

Instructor Demo of sodium acetate super-saturation

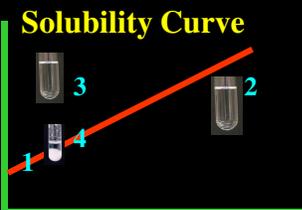

- 1 = Solid visible → Solution is saturated
- 2 = After heating: All solid dissolved → Solution unsaturated
- 3 = After slowly cooling: No solid visible → Supersaturated
- 4 = After Seeding: Solid visible → Saturated

Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

Sodium Acetate Demo Photos



Ambient

White Solid
Clear Liquid



~ 100 °C

No Solid
Clear Liquid



After Cooling

No Solid
Clear Liquid



After Seeding

White Solid
Clear Liquid

Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

Solubility of NaCl & KNO₃



NaCl

KNO₃



Ambient



~ 100 °C



After
Cooling



Ambient



~ 100 °C



After
Cooling

Large or Small
Increase or Decrease in solubility
Less visible “stuff” → greater the solubility


Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

Isolate 1.00 g of NaCl From 2 M Solution



- Calculate volume needed (See calculation slide)
- Pour 2 M solution into evaporating dish
- Cover with watch glass (limits spattering)
- Remove water by heating with a Bunsen Burner
- Weigh remaining solid

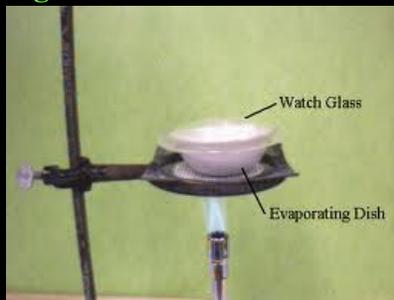

Hand On Set Up

Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

Volume of 2 M NaCl Solution Needed to Isolate 1.000 g NaCl

Determine Molecular Mass of NaCl

Na = 22.99

Cl = 35.45

58.44 g / mole

Dimensional analysis to solve for volume

$$1.000 \text{ g NaCl} \times \frac{1 \text{ mole NaCl}}{58.44 \text{ g}} \times \frac{1000 \text{ mL}}{2.000 \text{ mol}} = 8.560 \text{ mL}$$

The M means Moles per Liter or Moles per 1000 mL

Whenever you see M (Molarity), think moles / Liter or moles / 1000 mL

Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

Calculations

Mass of NaCl Solution:

Mass of evaporating dish , watch glass and NaCl Solution:
- Mass of evaporating dish and watch glass:
Mass of liquid:

Mass of NaCl Isolated:

Mass of evaporating dish , watch glass and NaCl Solid:
- Mass of evaporating dish and watch glass:
Mass of solid:

Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

Concentrations

% by Mass

$$\% \text{ (by mass)} = \frac{\text{grams solute}}{\text{grams solution}} \times 100$$



Molarity

$$\text{Molarity (M)} = \frac{\text{moles solute}}{\text{Liters solution}}$$



Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

Yields

$$\% \text{ Yield} = \frac{\text{Actual (Obtained in Experiment)}}{\text{Theoretical (Calculated Yield based on Stoichiometry)}} \times 100$$

$$\% \text{ Error} = \frac{\text{Actual Yield (g)} - \text{Theoretical Yield (g)}}{\text{Theoretical Yield (g)}} \times 100$$

% Error should be small and negative

Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

Conclusion



Describe the solubility of NaCl when the temperature changes

Delete incorrect term

Compare the solubility change for KNO₃ to the NaCl

Describe % yield

Describe % error



Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

Let's Boldly Go Explore Today's Lab



Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT