

% Water in Magnesium Sulfate

$\text{MgSO}_4 \cdot 7 \text{H}_2\text{O}$ → $\text{MgSO}_4 + 7 \text{H}_2\text{O}$

Copyright Larry P. Taylor, Ph.D. All Rights Reserved LPT

Hydrates

Absorb water from atmosphere
Water becomes associated with structure

$\text{MgSO}_4 \cdot 7 \text{H}_2\text{O}$

Copyright Larry P. Taylor, Ph.D. All Rights Reserved LPT

Hydrates

Compound + water → hydrate
reversible
Hydrate → compound + water

$\text{CuSO}_4 \cdot 5 \text{H}_2\text{O} \rightleftharpoons \text{CuSO}_4 + 5 \text{H}_2\text{O}$

$\text{MgSO}_4 \cdot 7 \text{H}_2\text{O} \rightleftharpoons \text{MgSO}_4 + 7 \text{H}_2\text{O}$

**Reactants and Products are chemically different:
Color change indicates chemical change
These reactions represent chemical changes**

Copyright Larry P. Taylor, Ph.D. All Rights Reserved LPT

Methane Hydrates (Clathrates)



**At cold temperatures:
Methane trapped by ice
Abundant in tundra and ocean bottoms
Global warming releasing the methane**







Copyright Larry P. Taylor, Ph.D. All Rights Reserved LPT

Desiccants

Compounds that absorb water to form hydrates









Used to protect variety of commercial products

Keep desiccants in containers until contents consumed


Copyright Larry P. Taylor, Ph.D. All Rights Reserved LPT



Desiccants

Five Common Types

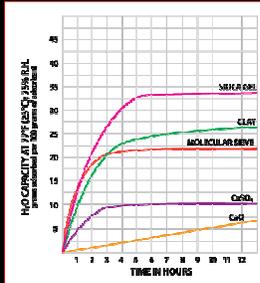

Time (h)	Molecular Sieves	Clay	Molecular Sieves	CaSO ₄	CaCl ₂
1	10	5	5	5	5
2	25	10	10	10	10
3	35	15	15	15	15
4	40	20	20	20	20
5	45	25	25	25	25
6	48	27	27	27	27
7	50	28	28	28	28
8	50	28	28	28	28
9	50	28	28	28	28
10	50	28	28	28	28
11	50	28	28	28	28
12	50	28	28	28	28

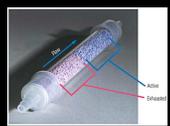

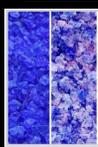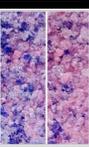Color Change shows absorption

Copyright Larry P. Taylor, Ph.D. All Rights Reserved LPT

Lab Desiccators

Used to Keep Sensitive Chemicals Dry






The Desiccator





Often stored under vacuum and sometimes in the cold (-78 °C)

Copyright Larry P. Taylor, Ph.D. All Rights Reserved LPT

Color of Many Gems From Hydrates





Never store in dehydrating conditions
Or
Underwater




Copyright Larry P. Taylor, Ph.D. All Rights Reserved LPT

Composition Calculations

Find the percent water present in the hydrate $\text{CuSO}_4 \cdot 5 \text{H}_2\text{O}$

Cu = 63.55		
S = 32.07	10 H = 10.08	
4 O = 64.00	5 O = 80.00	
Total = 159.62	Total = 90.08	
	Total = 249.70	

$\% \text{ Water} = \frac{90.08}{159.62 + 90.08} \times 100 = 36.08$
(249.70)

36.08 % = 0.3608



Copyright Larry P. Taylor, Ph.D. All Rights Reserved LPT

Composition Calculations

Find the percent water present in the hydrate $\text{MgSO}_4 \cdot 7 \text{H}_2\text{O}$

$$\text{Mg} = 24.31$$

$$\text{S} = 32.07$$

$$4 \text{ O} = 64.00$$

$$\text{Total} = 120.38$$

$$14 \text{ H} = 14.11$$

$$7 \text{ O} = 112.00$$

$$\text{Total} = 126.11$$



$$\% \text{ Water} = \frac{126.11}{120.38 + 126.11} \times 100 = 51.24$$

(246.49)





$$51.24 \% = 0.5124$$

Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

Naming Hydrates

Anhydrous (without water) name “• n H₂O’s”

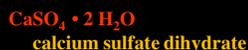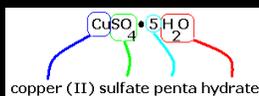

- Indicates distinct chemical entities held together

Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT


Desiccants keep my wraps dry



Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

% Water in Magnesium Sulfate Heptahydrate

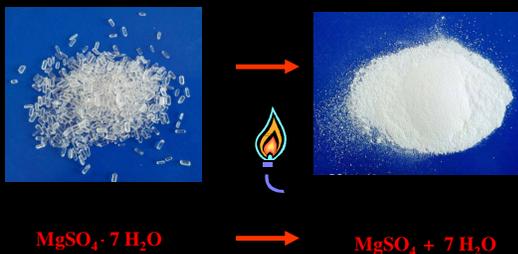

Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

Hydrate Lab

Purpose

Determine the percentage of water in a given hydrate

Procedure



Weigh materials "by difference"
Weight of evaporating dish, watch glass, & hydrate
- Weight of evaporating dish & watch glass

Weight of hydrate



Water driven away by heat
Watch glass minimizes splattering
Heat until all water is gone



Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

Calculations

Show all your work

Mass of a any substance (weighing by difference)

(Substance g + Container g) - Container g = Substance g

Theoretical Water loss: initial heptahydrate x % H_2O

Water Lost: Initial - final weight of the magnesium hydrate

Experimental % Water: $\frac{\text{mass H}_2\text{O lost}}{\text{mass initial heptahydrate}} \times 100$

Results

Tabulate the answers to your calculations

Conclusion

State % water in $\text{MgSO}_4 \cdot 7 \text{H}_2\text{O}$

Compare your experimental value to the theoretical

Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

Determine n



Calculate the value of n for $\text{MgSO}_4 \cdot n \text{H}_2\text{O}$

N is the ratio of moles water to moles anhydrous salt

Experiment measures grams ... need moles for this ratio

Convert grams water lost to moles (via molar mass of one H_2O)
Convert grams MgSO_4 remaining to moles (via molar mass MgSO_4)

$$n = \frac{\text{Moles water lost}}{\text{Moles anhydrous magnesium heptahydrate}}$$

n is closest small, whole number

Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT

Let's Boldly Go Explore Today's Lab



Copyright Larry P. Taylor, Ph.D. All Rights Reserved

LPT
